



Deleting and Adding a Probe

Deleting a Probe

1. Press Menu
2. Press Probe Settings
3. Change delete probe from off to on
4. Press List
5. On the list screen verify the green light flashing in the upper right corner (indicates Configuration file (CFU) is being sent to the ICC server).
6. Once the green light stops flashing, press the probe you would like to delete.
7. Verify you have chosen the correct probe and press the red delete button.
8. The screen will display “Delete Probe?” With the name of the probe to be deleted, if the correct probe has been chosen press delete again.
9. Press Menu
10. Press Probe settings
11. Change delete probe from on to off.
12. Press List
13. On the list screen verify the green light is flashing in the upper right corner.
14. Verify green light stops flashing, and the probe that you deleted is no longer showing up on the temperature list. (Should also check Temp Minder Online Probes page to confirm probe has been deleted).

Adding a Probe

1. Press Menu
2. Press Probe Settings
3. Change Probe Search from off to on.
4. Press List
5. On the list screen verify the green light flashing in the upper right corner
6. Once the green light stops flashing, pull the battery tab out of the new probe
7. Verify the screen displays “New Probe Found” and note the serial number of the probe you are adding appears on the screen.
8. Choose the position you would like to add the probe to by pushing the left and right arrows on the touch screen. (Note: If you accidentally choose to add the new probe to a position with an active probe the active probe will be replaced).
9. Press Menu
10. Press Probe Settings
11. Change probe search from on to off.
12. Press List
13. On the list screen verify the green light is flashing in the upper right corner.
14. Verify the green light stops flashing, and that the probe you added is on the temperature list. (Should also check Temp Minder Online Probes page to confirm new probe is added).
15. Hang the probe in Fridge Or Freezer

